

NCL (New Cereal Leaf-Prescott Liquid)

Protozoa

Medium	per litre
Cerophyll, cereal grass leaves *	1.0 g
Prescott's & James's Solution (PJ - see recipe below)	1.0 litre

Bring PJ to the boil then add cereal leaves. Continue to boil for 5 minutes. Allow to cool then restore volume to 1 litre with deionized water. Filter through GF/C paper and autoclave at 15 psi for 15 minutes.

Supply

* Ward's Natural Science, PO Box 92912, Rochester, NY 14692-9012, USA
<http://wardsci.com>

PJ (Prescott's & James's Solution)

Stocks	per 100 ml
(1) CaCl ₂ .2H ₂ O	0.43 g
KCl	0.16 g
(2) K ₂ HPO ₄	0.51 g
(3) MgSO ₄ .7H ₂ O	0.28 g

Medium	per litre
Stock solutions 1 - 3	1.0 ml each

Make up to 1 litre with deionized water. Autoclave at 15 psi for 15 minutes.