

ASW 225 + Barley (Artificial seawater with salinity of 225 psu)

Stocks	per 1000 cm ³
(1) NaCl	212.6 g
KCl	3.8 g
MgCl ₂ .6H ₂ O	13.45 g
MgSO ₄ .7H ₂ O	1.65 g
CaCl ₂ .2H ₂ O	0.65 g

Medium

Add 20 ml of the autoclaved artificial seawater with 255 psu (stock 1) and 4 autoclaved barley grains to a T-25 tissue culture flask (Corning 25100).

Used for CCAP 1675/1 *Trimyema koreanum*