

MASM (Modified Artificial Seawater Medium)

Stocks

per 100 ml

(1) MgSO ₄ .7H ₂ O	24.4 g
(2) KCl	6.0 g
(3) NaNO ₃	10.0 g
(4) CaCl ₂ .2H ₂ O	3.0 g
(5) KH ₂ PO ₄	0.5 g
(6) NH ₄ Cl	2.67 g
(7) Trace Elements (PIV):	
Ensure elements are added in the following sequence:	
Na ₂ EDTA	0.75 g
FeCl ₃ .6H ₂ O	0.097 g
MnCl ₂ .4H ₂ O	0.041 g
ZnCl ₂ .6H ₂ O	0.005 g
CoCl ₂ .6H ₂ O	0.002 g
Na ₂ MoO ₄ .2H ₂ O	0.004 g

Once elements are dissolved autoclave at 15 psi for 15 minutes.

Per 100 ml

- | | |
|---|--------|
| (8) Vitamin B ₁ (Thiamine hydrochloride) | 0.12 g |
| Filter sterile | |
| (9) Vitamin B ₁₂ (Cyanocobalamin) | 0.1 g |
| Take 1 ml of this solution and add 99 ml Deionised water. Filter sterile. | |
| (10) Soil extract 2 – SE2 (see overleaf) | |

Medium

per litre

Tris	1.0 g
NaCl *	30 g
Stock solution 1 - 5	10 ml each
Stock solution 6	1 ml
Stock solution 7 (Trace element)	6 ml
Stock solutions 8 - 9	1 ml each
Stock solution 10 (SE2)	30 ml

Make up to 1 litre with distilled water and adjust to pH **8.0** with 1M NaOH or 1M HCl.
Autoclave at 15 psi for 15 minutes.

* For brackish organisms, take 15 g of NaCl instead of 30 g (BW/MASM)

Reviewed: 10th August 2020

Created on: 05 Nov 2019	CCAP (Culture Collection of Algae and Protozoa), SAMS Ltd, Scottish Marine Institute, Oban, Argyll, PA37 1QA, UK Tel: +44 (0)1631 559000 Fax: +44 (0)1631 559001 Email: ccap@sams.ac.uk Web: www.ccap.ac.uk	2 Pages
----------------------------	--	---------

SE2 (Soil Extract 2)

Freshwater and terrestrial protozoa

Preparing the soil

Site selection for a good soil is very important and for most purposes a soil from undisturbed deciduous woodland is best. Sites to avoid are those showing obvious signs of man's activity and particular care should be taken to avoid areas where fertilizers, crop sprays or other toxic chemicals may have been used.

A rich loam with good crumb structure should be sought. Stones, roots and larger invertebrates should be removed during an initial sieving through a 1 cm mesh. The sieved soil should be spread to air dry and handpicked for smaller invertebrates and roots. It should be turned periodically and picked over again. When dry sieve through a finer mesh (2-4 mm) and store in an airtight container away from light and heat.

Medium

Soil is prepared as above. Air-dried soil and twice its volume of supernatant distilled water are autoclaved together at 15 psi for 2 hours and left to cool. The supernatant is then decanted and then distributed to containers in volumes suitable for making up batches of media. The aliquots and their containers are autoclaved for an appropriate length of time (e.g. 1 litre or less for 15 minutes) and are then kept in a cool place (e.g. a refrigerator) until required.

Reviewed: 10th August 2020

Created on: 05 Nov 2019	CCAP (Culture Collection of Algae and Protozoa), SAMS Ltd, Scottish Marine Institute, Oban, Argyll, PA37 1QA, UK Tel: +44 (0)1631 559000 Fax: +44 (0)1631 559001 Email: ccap@sams.ac.uk Web: www.ccap.ac.uk	2 Pages
----------------------------	---	---------