

MY75S (Malt & Yeast Extract - 75% Seawater Agar)

Marine Protozoa

Medium**per litre**

Natural Filtered seawater	750.0 ml
Deionized water	250.0 ml
Malt extract (Oxoid L39) *	0.1 g
Yeast extract (Oxoid L21) *	0.1 g
Bacteriological Agar (Oxoid L11) *	15.0 g

Make up the 75% seawater, heat, (do not bring to boil) then add and dissolve the malt extract and yeast extract using a magnetic stirrer. Cool slightly then add the agar and shake to disperse. Autoclave at 15 psi for 15 minutes.

Supply

* Unipath Ltd, Wade Road, Basingstoke, Hants., RG24 0PW, UK

Reviewed: 5th October 2020