

## PAS (Page's Amoeba Saline Solution)

Constituent of protozoa media

### Stocks

per 500 ml

(1) NaCl	12.0 g
MgSO <sub>4</sub> .7H <sub>2</sub> O	0.40 g
CaCl <sub>2</sub> .2H <sub>2</sub> O	0.60 g
(2) Na <sub>2</sub> HPO <sub>4</sub>	14.20 g
KH <sub>2</sub> PO <sub>4</sub>	13.60 g

### Medium

per litre

Stock solutions 1 & 2	5.0 ml each
-----------------------	-------------

Make up to 1 litre deionised water. Autoclave at 15 psi for 15 minutes.

Reviewed: 10<sup>th</sup> August 2020