

PJ (Prescott's & James's Solution)

Freshwater protozoa

Stocks**per 100 ml**

(1) CaCl ₂ .2H ₂ O	0.43 g
KCl	0.16 g
(2) K ₂ HPO ₄	0.51 g
(3) MgSO ₄ .7H ₂ O	0.28 g

Medium**per litre**

Stock solutions 1 - 3	1.0 ml each
-----------------------	-------------

Make up to 1 litre with deionized water. Autoclave at 15 psi for 15 minutes.

Reviewed: 10th August 2020