

PP (Proteose Peptone Medium)

Freshwater algae

Stocks

per litre

(1) K ₂ HPO ₄	1.0 g
(2) MgSO ₄ .7H ₂ O	1.0 g
(3) KNO ₃	10.0 g

Medium

per litre

Proteose peptone (Oxoid L85) *	1.0 g
Stock solutions 1 - 3	20.0 ml each

The stock solutions are those for SES. Make up to 1 litre with distilled water. For agar add 15g per litre Bacterial Agar. Autoclave at 15 psi for 15 minutes.

Supply

* Unipath Ltd, Wade Road, Basingstoke, Hants, RG24 0PW, UK

Reviewed: 10TH August 2020