

PPY (Proteose Peptone Yeast Extract Medium)

Freshwater protozoa

Medium

per litre

Proteose peptone (Oxoid L85) *	20.0 g
Yeast extract (Oxoid L21) *	2.5 g

Make up to 1 litre with deionized water. Autoclave at 15 psi for 15 minutes.

Note. For test tubes, only use paper or pink silicone bungs.

Supply

* Unipath Ltd, Wade Road, Basingstoke, Hants, RG24 0PW, UK

Reviewed: 10th August 2020